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Strumento: EASY-Spray Source Kit (NG)  
 ID di serie: ES041794

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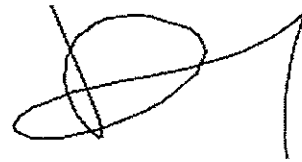
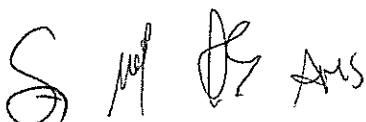
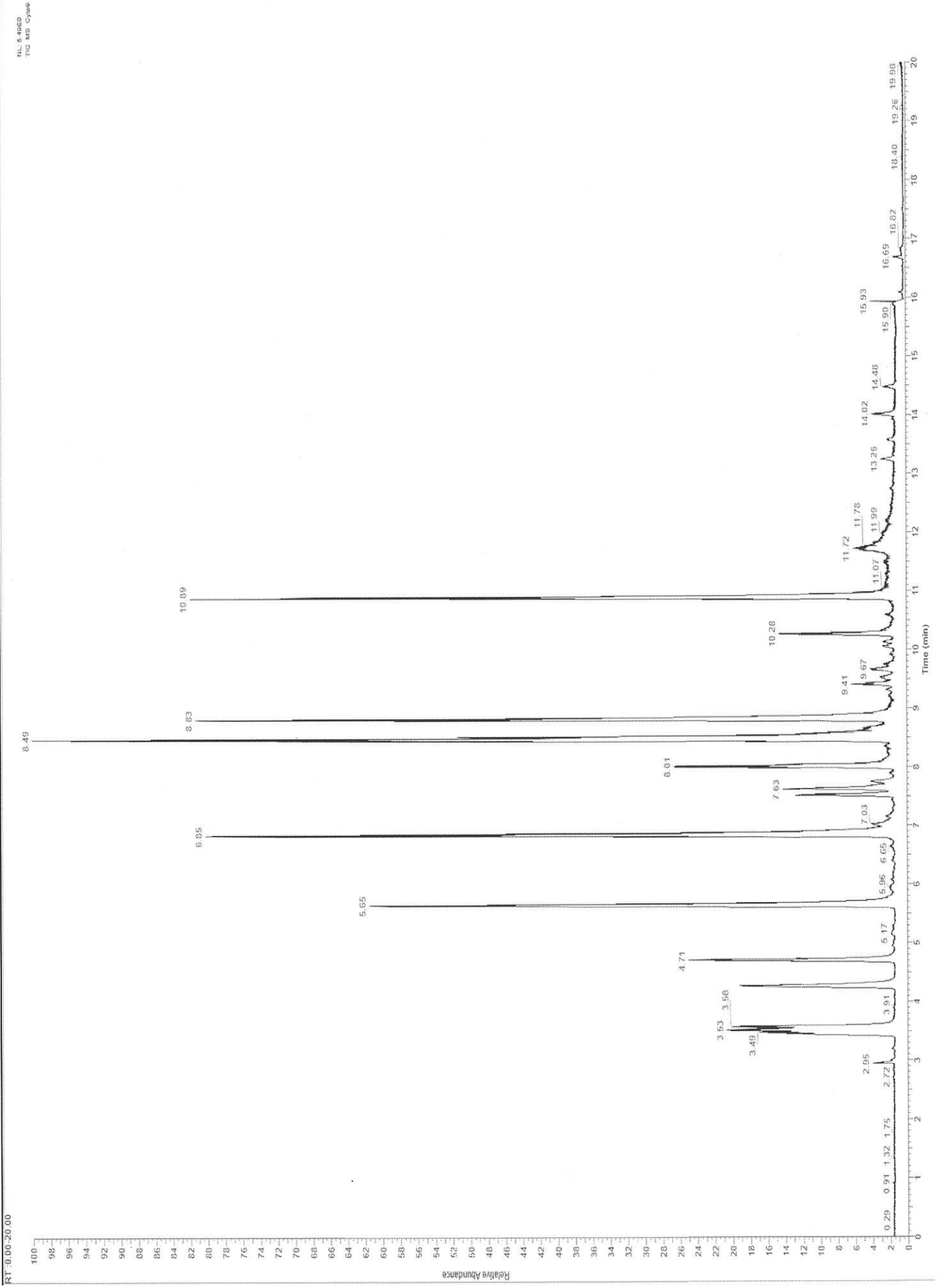
Data: 26-Mar-2024

MS: 6.4020  
Thu, Mar 07, 2006

*Handwritten signatures and initials in blue ink.*

# Diagnostic Report

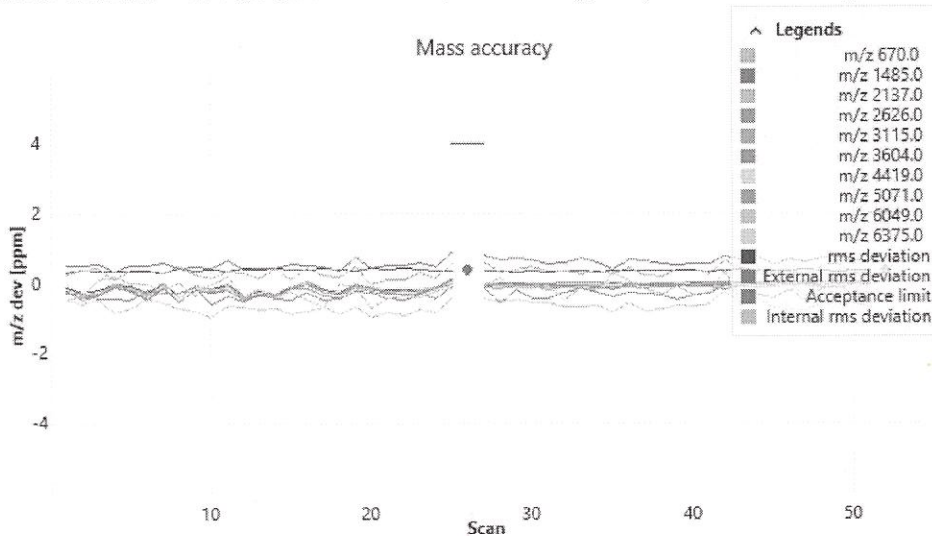
Date & Time Monday, March 25, 2024 04:39:20 PM  
Instrument Model Orbitrap Exploris 480  
Instrument Serial MA11076C  
Software Version 4.2.362.42

Name	Result	Comment
High m/z AHFP Spectral Mass Accuracy Calibration (pos)	Passed	External rms deviation = 0.38 ppm (without lock mass), acceptance limit = 4.00 ppm Internal rms deviation = 0.34 ppm (with lock mass m/z 1485.02082), acceptance limit = 2.00 ppm

## High m/z AHFP Spectral Mass Accuracy Calibration (pos)

External rms deviation (without lock mass): 0.38 ppm is in spec. Value does not exceed 4 ppm  
Internal rms deviation (with lock mass m/z 1485.02082): 0.34 ppm is in spec. Value does not exceed 2 ppm

High m/z AHFP Spectral Mass Accuracy Calibration (pos) passed:  
External rms deviation = 0.38 ppm (without lock mass), acceptance limit = 4.00 ppm  
Internal rms deviation = 0.34 ppm (with lock mass m/z 1485.02082), acceptance limit = 2.00 ppm


## System Configuration

Name	Value
Ion Source Type	ESI
Internal Calibration (EASY-IC) Source	enabled
Options	EASY-IC, BioPharma

## Embedded System Information

Label	Version
Instrument Embedded Software	4.2.319.30-SP4
Instrument Hardware Hash	8e3c4ae04ad27dab2fae82686fbb05a8830272f8

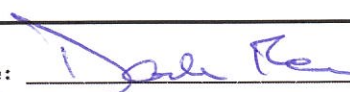Serial Number: MA11076C

Signature: 

Monday, March 25, 2024 04:40:22 PM

*So M AS AM*

# Diagnostic Report

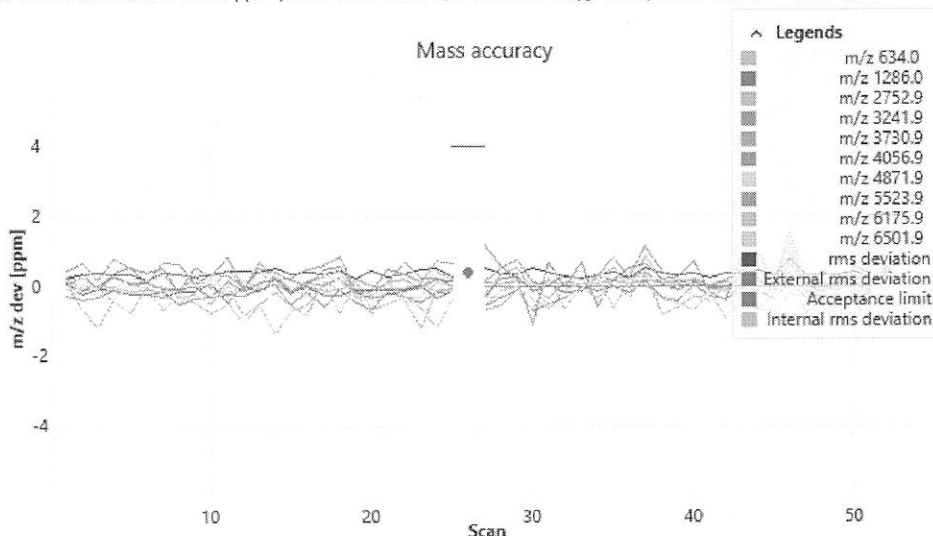
Date & Time      Monday, March 25, 2024 04:45:16 PM  
 Instrument Model    Orbitrap Exploris 480  
 Instrument Serial    MA11076C  
 Software Version    4.2.362.42

Name	Result	Comment
High m/z AHFP Spectral Mass Accuracy Calibration (neg)	Passed	External rms deviation = 0.37 ppm (without lock mass), acceptance limit = 4.00 ppm Internal rms deviation = 0.38 ppm (with lock mass m/z 1285.95461), acceptance limit = 2.00 ppm

### High m/z AHFP Spectral Mass Accuracy Calibration (neg)

External rms deviation (without lock mass): 0.37 ppm is in spec. Value does not exceed 4 ppm  
 Internal rms deviation (with lock mass m/z 1285.95461): 0.38 ppm is in spec. Value does not exceed 2 ppm

High m/z AHFP Spectral Mass Accuracy Calibration (neg) passed:  
 External rms deviation = 0.37 ppm (without lock mass), acceptance limit = 4.00 ppm  
 Internal rms deviation = 0.38 ppm (with lock mass m/z 1285.95461), acceptance limit = 2.00 ppm

### System Configuration

Name	Value
Ion Source Type	ESI
Internal Calibration (EASY-IC) Source	enabled
Options	EASY-IC, BioPharma

### Embedded System Information

Label	Version
Instrument Embedded Software	4.2.319.30-SP4
Instrument Hardware Hash	8e3c4ae04ad27dab2fae82686fbb05a8830272f8

Serial Number: MA11076C

Signature:   
 Monday, March 25, 2024 04:47:20 PM

SA    MM    AS    JMS